

CERTIFICATO
ILVO

AnticFast® 3 in 1


Test rapidi certificati ILVO per ricerca di

Cat.# MZLFBTC Beta-lattamici+Tetracicline+Cefalexina

Cat.# MZLFBTS Beta-lattamici+Tetracicline+Sulfamidici




PROCEDURA

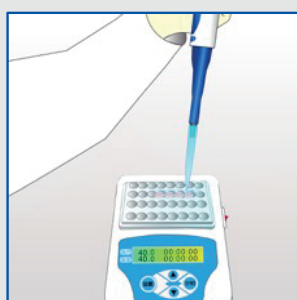


Aggiungere 200 µl di latte



Incubare per il tempo indicato in tabella

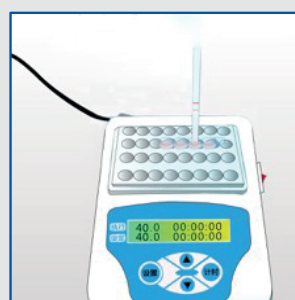

Sviluppare per 5 minuti a 40 C°

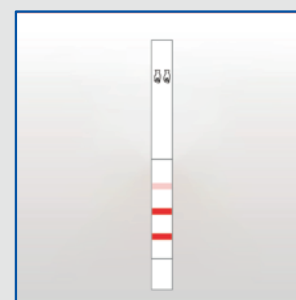

Interpretazione visiva


INTERPRETAZIONE VISIVA DEI RISULTATI

Cat.# MZLFBTC

T1: Tetracicline	3 minuti
T2: Beta-lattamici	
T3: Cefalexina	

Cat.# MZLFBTS

T1: Sulfamidici	2 minuti
T2: Tetracicline	
T3: Beta-lattamici	

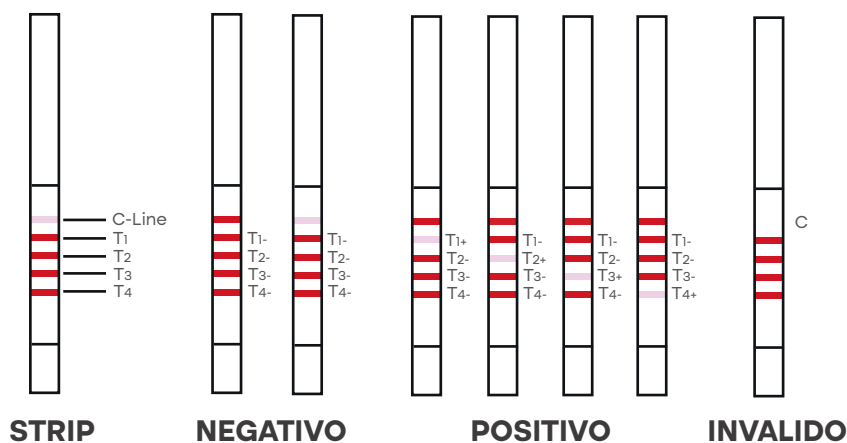



MATERIALE CONTENUTO NEL KIT

- Manuale
- 96 tests/kit, 12 tubi da 8 test LFD
- 96 pozzetti
- Controllo Positivo/negativo
- Pipetta
- Puntali

Scadenza	Conservazione
12 mesi	2-8 °C

LIMITI DI DETECTION

Penicilline	LOD (ppb)	Tetracicline (T)	LOD (ppb)	Sulfamidici (S)*	LOD (ppb)
Amoxicillin	< EU MRL (µl/Kg)	Chlortetracycline	< EU MRL (µl/Kg)	Sulfamethoxyprizinum	< EU MRL (µl/Kg)
Ampicillin	< EU MRL (µl/Kg)	Doxycycline	< 18	Sulfametoxydiazine	< EU MRL (µl/Kg)
Benzathine	< 5	Oxytetracycline	< EU MRL (µl/Kg)	Sulfamonomethoxine	< EU MRL (µl/Kg)
Cloxacillin	< EU MRL (µl/Kg)	Tetracycline	< EU MRL (µl/Kg)	Sulfaquinoxaline	< EU MRL (µl/Kg)
Dicloxacillin	< EU MRL (µl/Kg)	Sulfamidici (S)*	< EU MRL (µl/Kg)	Sulfisomidine	< EU MRL (µl/Kg)
Nafcillin	< EU MRL (µl/Kg)	Sulfabenzamide	< EU MRL (µl/Kg)	Sulfisoxazole	< EU MRL (µl/Kg)
Oxacillin	< EU MRL (µl/Kg)	Sulfachlorpyridazine	< EU MRL (µl/Kg)		
Penicillin G	< EU MRL (µl/Kg)	Sulfaclozine	< EU MRL (µl/Kg)		
Cefalosporine	LOD (ppb)	Sulfadiazine	< EU MRL (µl/Kg)		
Cefquinome	< EU MRL (µl/Kg)	Sulfadimethoxine	< EU MRL (µl/Kg)		
Ceftiofur	< EU MRL (µl/Kg)	Sulfadoxine	< EU MRL (µl/Kg)		
Cefalonium	< EU MRL (µl/Kg)	Sulfaethoxypridazine	< EU MRL (µl/Kg)		
Cephapirin	< EU MRL (µl/Kg)	Sulfamerazine	< EU MRL (µl/Kg)		
Cefacetrile	< EU MRL (µl/Kg)	Sulfamethazine	< EU MRL (µl/Kg)		
Cefazolin	< EU MRL (µl/Kg)	Sulfamethazole	< EU MRL (µl/Kg)		
Cefoperazone	< EU MRL (µl/Kg)	Sulfamethizole	< EU MRL (µl/Kg)		
Cefalotin*	< 20	Sulfamethoxazole	< EU MRL (µl/Kg)		
Cefalexina**	< EU MRL (µl/Kg)	Sulfamethoxypridazine	< EU MRL (µl/Kg)		

*solo nel kit MZLFBTS

**solo nel kit MZLFBTC